## **FIERCE LA-ICP-MS sample holders:**

- Slide sample holder (ideal for thin sections, polished blocks, round and rectancular)
- Round sample holder (ideal for standard sized epoxy mounts)
- Universal sample holder (ideal for oddly shaped samples, even >100mm)

**Principle**: all samples and standards are pushed upwards relative to a reference surface with springs. In order to guarantee a flat and smooth surface (which improves visibility and data quality) polishing of samples is advised.







## Universal sample holder

This sample holder is perfectly suited to accomodate unregularly shaped samples



Samples are mounted and held in place with clamps and spacers. Small springs beneath the sample push the sample upwards.